

Install PSI Driver

1. Plug the USB connector to the PC and then switch the power on.



2. Found New Hardware Wizard appears.



Click Next button.

3.




Select Search for a suitable driver for my device (recommended) and click Next button.

4.



Click Browse... button and locate the directory where the PSI USB Driver files are stored. For example: C:\Install\PSI\FluorCam\PSI\USB_Driver\V1.0.0.3. Be sure that you have version 1.0.0.3 of PSI USB Driver.

5.

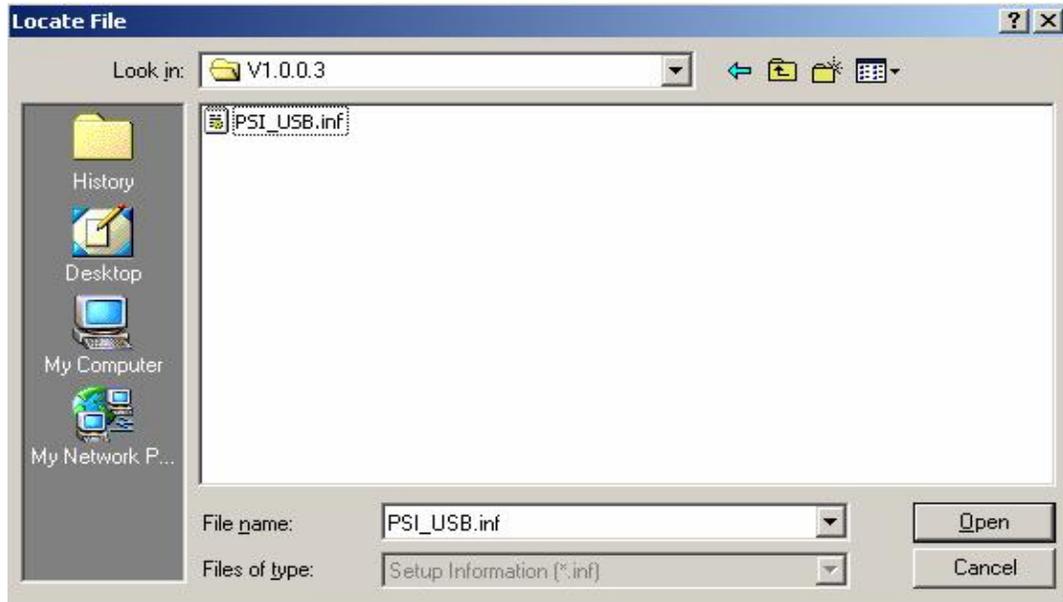


Find PSI_USB.inf file and click Open button.

6.



Click OK button.

7.

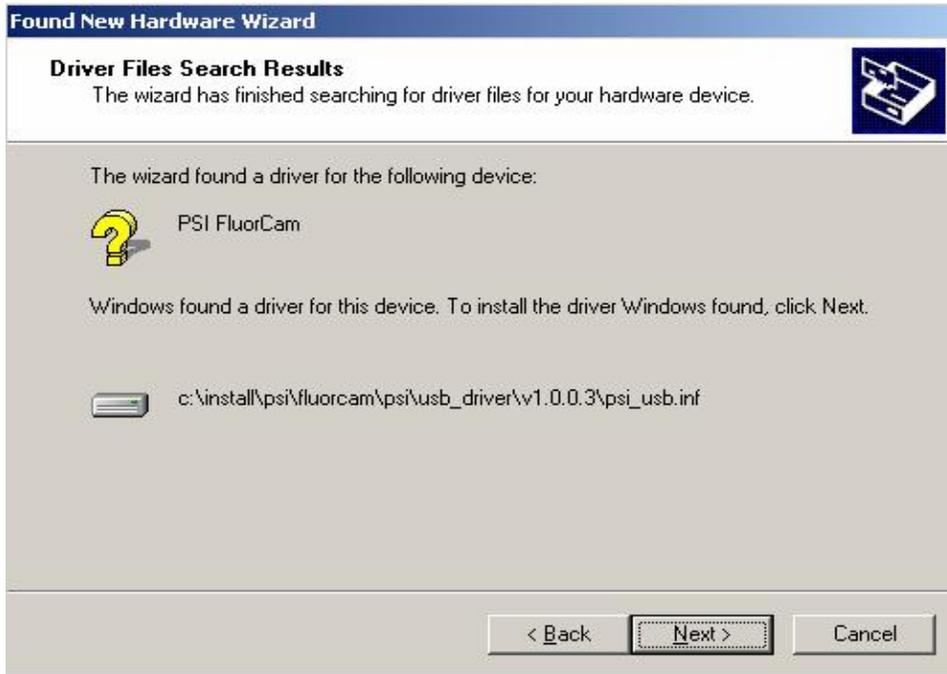


Click Next button.

8.




Click Finish button.

9. Right click on My Computer and select Properties. System Properties Window appears.

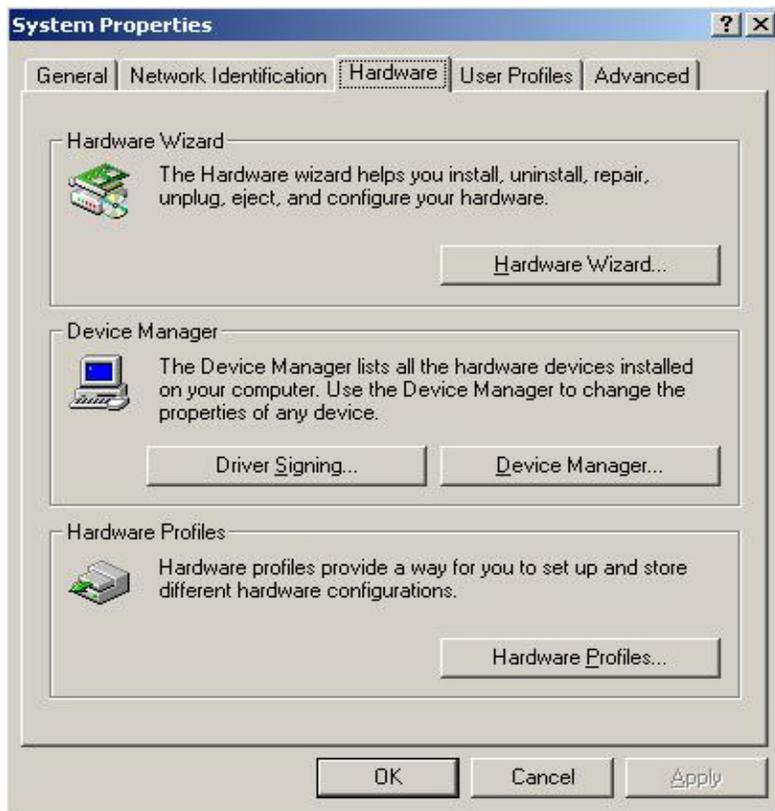


Go to Hardware page and click Device Manager button.

10.

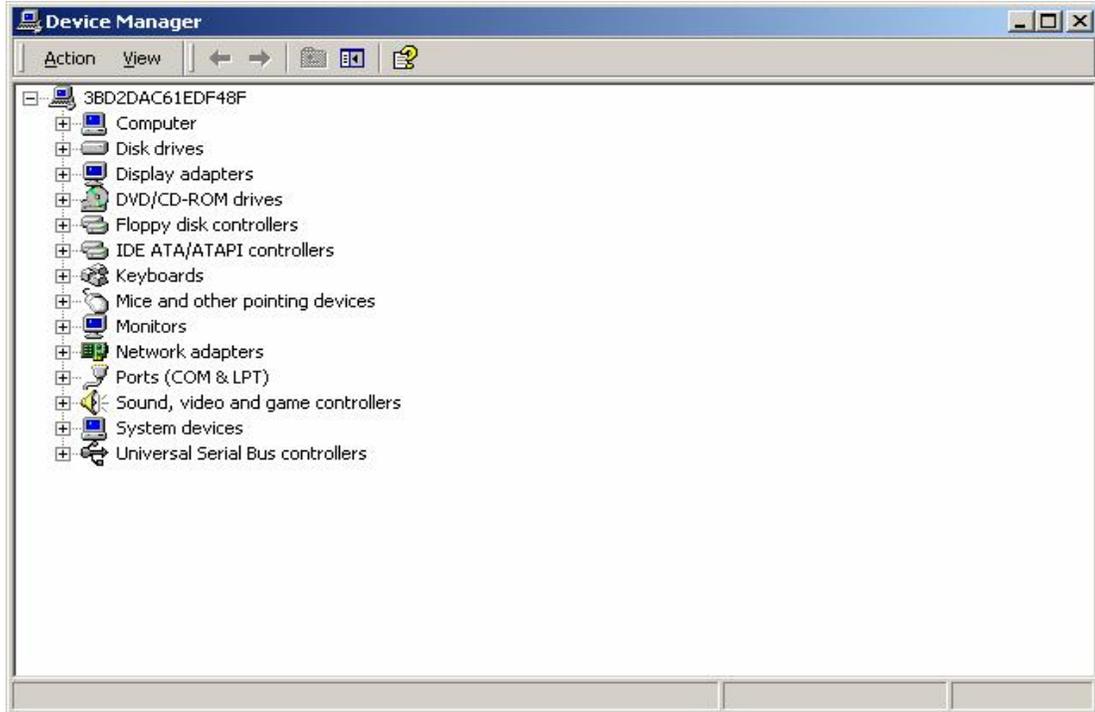

Click + in the Universal Serial Bus controllers entry.

11.

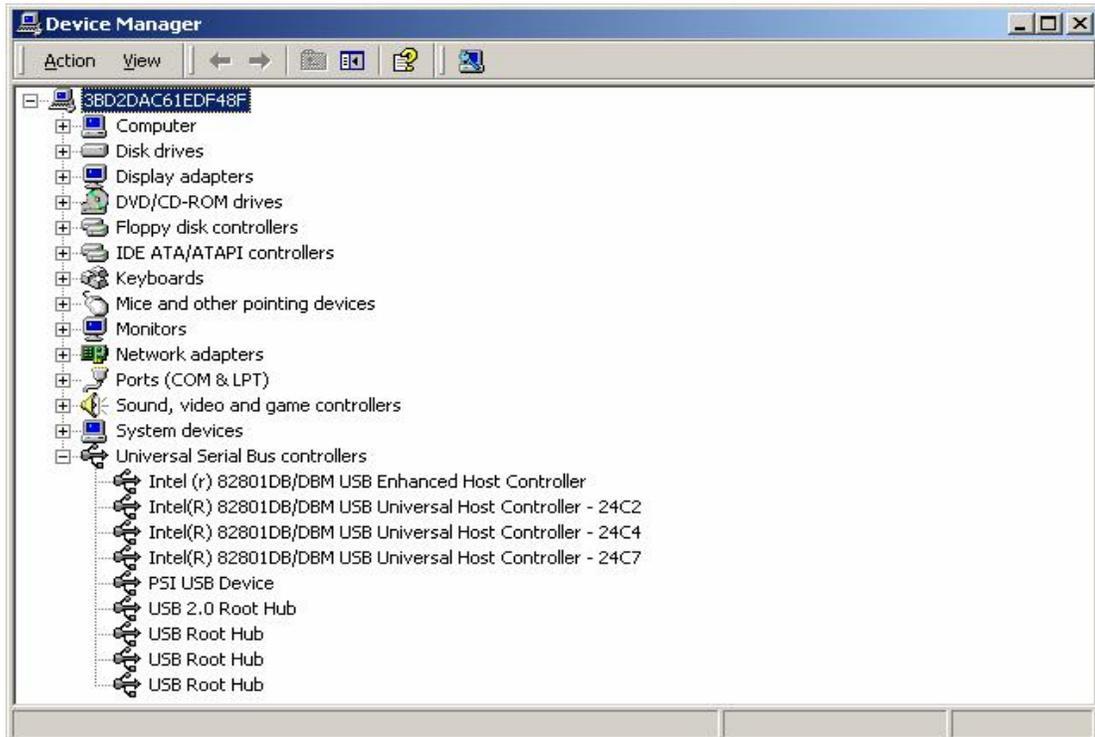

You have to see PSI USB Device entry without yellow exclamation mark.
In this case the PSI USB Device is installed correctly.